# Time-Resolved Radioluminescence of Perovskite Scintillators

RH\_P15; Tommy Loan

### Introduction

Scintillators are materials that convert X-rays into UV or visible radiation for detection with a photodetector or imaging sensor. Such materials are in high demand for a diverse range of applications, including medical imaging, security checks, and scientific research. Metal halide perovskites such as CsPbBr<sub>3</sub> nanocrystals (NCs) have attracted attention due to their bright radioluminescence (RL), high X-ray absorbance, and low detection limit.

This Research Highlight presents a study published in 2023 in ACS Applied Nano Materials by researchers at the University of Surrey.<sup>1</sup> In this work, the scintillation properties of CsPbBr<sub>3</sub> NCs were characterised with an Edinburgh Instruments FLS1000 Photoluminescence (PL) Spectrometer equipped with the XS1 RL chamber

## Materials and Methods

Batches of CsPbBr<sub>3</sub> NCs were synthesised with different volumes (0.0, 0.05, 0.1, 0.2, 0.4, 0.8 mL) of an oleylamine (OAM) surfactant and were named accordingly (OAM 0.0, OAM 0.4 etc.). NCs were then loaded into poly(methyl methacrylate) (PMMA) plastic discs at a range of concentrations (1, 2, 4, and 10%). RL emission of the CsPbBr<sub>3</sub> NCs / PMMA discs was measured with an FLS1000 spectrometer and XS1 RL chamber (Figure 1). Steady-state RL emissions of the samples were measured under excitation by a 60 kV X-ray with a tube current of 100  $\mu$ A. Time-resolved RL decays were acquired using TCSPC with a laser-excited 40 kV pulsed X-ray source.



Figure 1 Edinburgh Instruments XS1 RL chamber.

### Results

The scintillation properties of the discs were characterised in the XS1 and are shown in Figure 2. A 4% OAM 0.4-loaded PMIMA disc (PMIMA OAM 0.4) produced an emission band at 536 nm (Figure 2a). It was found that RL intensity of the discs increased with both OAM concentration (Figure 2b) and NC loading (Figure 2c).

EDINBURGH INSTRUMENTS

The RL decay of the 4% CsPbBr<sub>3</sub> NC/PMMA disc was measured using a pulsed X-ray source, revealing a triexponential decay with an average RL lifetime of 29.4 ns (Figure 2d), which is significantly faster than commercial scintillators such as CsI:TI (~680 ns decay time).

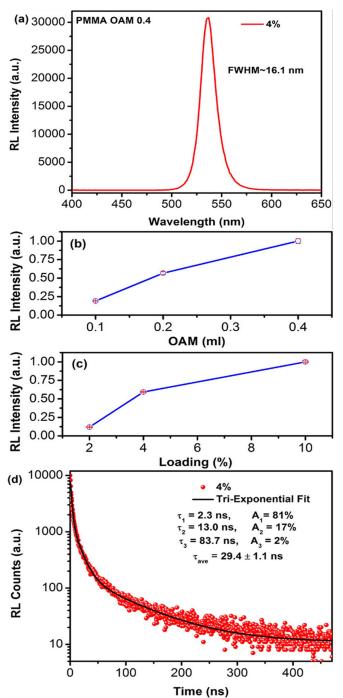


Figure 2 (a) RL emission spectrum of PMMA disc loaded with 4% OAM 0.4. (b) Plot of RL emission against OAM volume for CsPbBr<sub>3</sub> NC/PMMA disc. (c) Plot of RL emission against NC loading of PMMA discs. (d) RL lifetime decay and fitting of a 4% CsPbBr<sub>3</sub> NC/PMMA disc. The image is reprinted from Ghosh et al.<sup>1</sup>, Copyright (2024), with permission from American Chemical Society.

#### **RESEARCH HIGHLIGHT**

# Time-Resolved Radioluminescence of Perovskite Scintillators

RH\_P15; Tommy Loan



## Conclusions

In this Research Highlight, it was shown how the FLS1000 equipped with the XS1 RL chamber can be used to characterise novel scintillator materials. The intensity of RL emission of the scintillators could be tuned with the surfactant concentration and NC loading. The materials exhibited a maximum RL emission band at 536 nm with a rapid lifetime of 29.4 ns, demonstrating its ability to act as an X-ray detector.

#### References

 Ghosh, J. et al. Surfactant-Dependent Bulk Scale Mechanochemical Synthesis of CsPbBr<sub>3</sub> Nanocrystals for Plastic Scintillator-Based X-ray Imaging. ACS Appl. Nano. Mater. 6, (2023).

### **Full Publication**

The full article was published in ACS Applied Nano Materials. It is available at: https://pubs.acs.org/doi/full/10.1021/acsanm.3c02531.



For more information, please contact:

+44 (0) 1506 425 300 sales@edinst.com www.edinst.com